



YamayBio

Tris/Glycine/SDS Buffer Instant Granules

USER GUIDE

For research use only.
Not for use in diagnostic procedures.



Contents and Storage

Tris/Glycine/SDS Buffer Instant Granules

MS8131

10 sticks (1 L×10)

Storage: Store at room temperature for 36 months. The product is shipped at room temperature.

Introduction

This product is supplied as white or pale yellow instant granules composed of Tris, glycine, and SDS. Each stick is sufficient to prepare 1 L of 1× Tris/glycine/SDS electrophoresis buffer, with final concentrations 25 mM Tris, 192 mM glycine, and 0.1% SDS.

Quick Start Protocol

Dissolve one stick in distilled water and bring the final volume to **1 L** to prepare 1× Tris/glycine/SDS electrophoresis buffer.

