



YamayBio

Transfer Buffer Instant Granules

USER MANUAL

For research use only.
Not for use in diagnostic procedures.

Contents and Storage

Transfer Buffer Instant Granules

MS8134

10 sticks (1.2 L×10)

Storage: Store at room temperature for 36 months. The product is shipped at room temperature.

Introduction

This product is supplied as white or pale yellow instant granules. The main components are Tris and glycine. Each stick prepares 1.2 L of 1× transfer buffer for wet transfer applications in Western blotting.

Quick Start Protocol

1. Add approximately **1 L** of **distilled water** to a beaker. Slowly add one stick of Transfer Buffer Instant Granules, stirring continuously until the granules are fully dissolved.
2. Add **200 mL of anhydrous ethanol or methanol** to the resulting solution. Mix thoroughly to obtain 1× wet transfer buffer.