



YamayBio

Tricine SDS Running Buffer Instant Granules

USER GUIDE

For research use only.
Not for use in diagnostic procedures.



Contents and Storage

Tricine SDS Running Buffer Instant Granules

MS8136 10 sticks (500mL×10)

Storage: Store at room temperature for 36 months. The product is shipped at room temperature.

Introduction

This product is supplied as white or pale yellow instant granules. Each stick is sufficient to prepare 500 mL of 1× Tricine SDS electrophoresis buffer.

When used for Tricine-SDS-PAGE, separate anode and cathode buffers are not required.

Quick Start Protocol

Dissolve one stick in distilled water and bring the final volume to **500 mL** to obtain 1× Tricine SDS electrophoresis buffer.

